




SPure™ HC Filters with Push-to-Connect Fittings

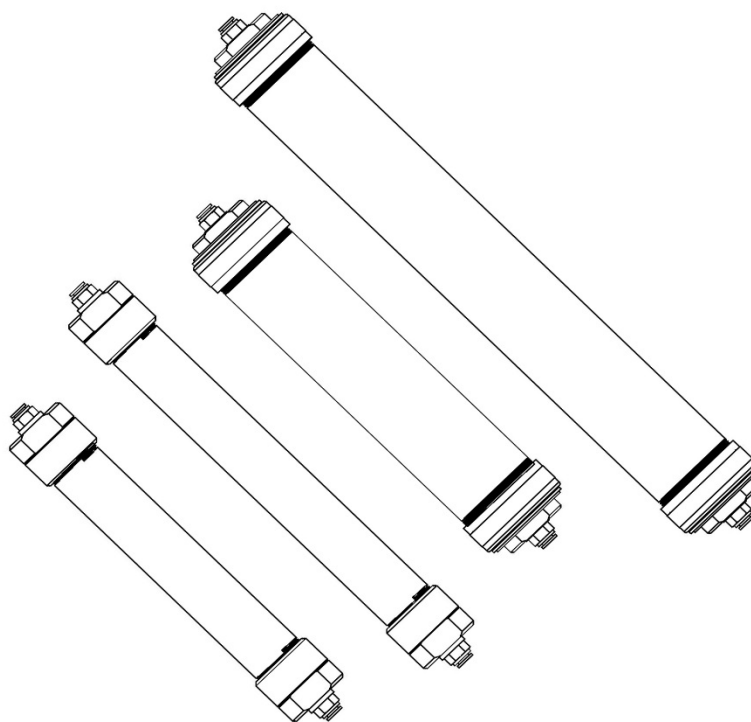



Warnings

Warnings must be read carefully and understood. Improper use of this product can cause harm or death to personnel and damage to property!



Wear eye protection and use caution when working with pressurized systems.

Maximum pressure rating for SPure HC Filters: 6.9 bar / 100 psi.


When using a high-pressure gas source overpressure protection must be provided.

Do not expose the protective outer plastic tube to solvents.

Not for use with hydrogen or hazardous gases.
Consult local regulations and your company's safety procedures.

Product Numbers

Size	For 1/4" Tubing	For 6mm Tubing	Maximum Pressure
88 cc	202326-4-PC	202326-6-PC	6.9 bar / 100 psi
116 cc	202326L-4-PC	202326L-6-PC	6.9 bar / 100 psi
240 cc	202326D-4-PC	202326D-6-PC	6.9 bar / 100 psi
400 cc	202326XL-4-PC	202326XL-6-PC	6.9 bar / 100 psi

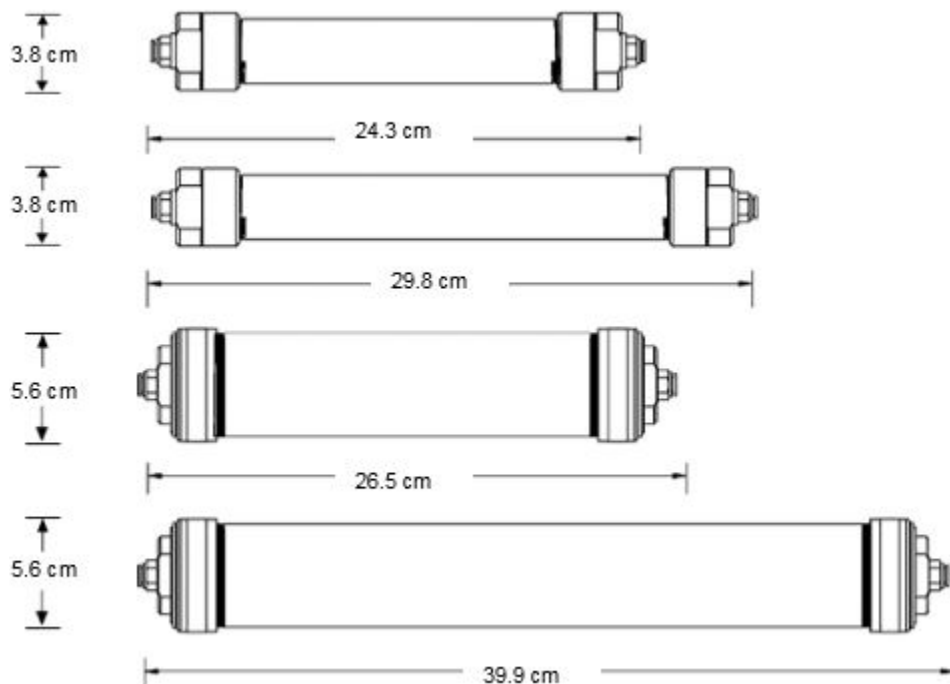

From top to bottom: 88 cc, 116 cc, 240 cc, 400 cc.

Filter Setup and Use

Installation for Push-to-Connect Fittings

1. Check the package contents for damage. Contact your supplier if any items are damaged or missing.
2. The SPure Filter is delivered in the active state, filled with helium. It is ready for use, but it is recommended that the trap be purged with the gas to be purified.
3. Make sure the ends of the plastic tubing are clean and free of scratches. If it is damaged, simply cut a small piece off the end.
4. Remove the shipping plug from the inlet end of the trap by depressing the locking tab and pulling on the plug.
5. Insert the tube into the fitting until it bottoms out. Make sure the tube is fully inserted.
6. Gently pull back on the tube to make sure the locking tab is engaged.
7. Repeat for the outlet end of the trap.
8. Purge the entire system until a total of 1 to 2 liters of gas has been flushed through. Maintain a low purge rate of 10 to 30 cc/min
9. Leak check all connections.
10. If leaks are present, remove the tube by pressing down on the locking tab, and simultaneously pulling up on the tube. Trim the end of the tube and re-insert.



Trajan Scientific Americas Inc
2601 Technology Drive,
Louisville, KY 40299, USA
Tel: 502 491 6300
Fax: 502 491 3390
usa@trajanscimed.com



Recycling

Please contact CRS for information regarding recycling and reconditioning gas purifier products.

© Trajan Scientific Americas Inc
09/2022 995250 _RevB