


## SPure™ H<sub>2</sub>O/HC filters

### Quick connect fittings

- **Color Disclaimer:** This product is intended to be yellow. Slight variations in shade may occur due to the manufacturing process and differences in individual color perception.

### Introduction

Pure gas is critical requirement in gas chromatography, spectroscopy, optics, lithography, and numerous other applications in manufacturing and analytical laboratories. The SPure™ H<sub>2</sub>O/HC filters removes moisture, oil, dust and hydrocarbons (C5 and heavier) from dry air and inert gases.

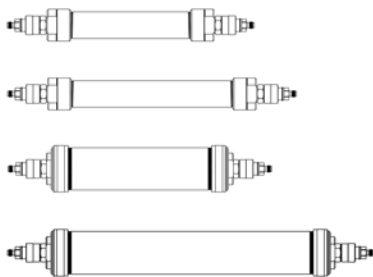

### Important



Warnings must be read carefully and understood. Improper use of this product can cause harm or death to personnel and damage to property!

- Wear eye protection and use caution when working with pressurized systems.
- Maximum pressure is 6.9 bar / 100 psi.
- When using a high-pressure gas source over-pressure protection must be provided.
- Do not use with supply gas H<sub>2</sub>O concentration greater than 0.1%. Contains molecular sieve which may become very hot in contact with liquid water. Do not use with liquids.
- Do not expose the protective outer plastic tube to solvents.
- Not for use with hydrogen or hazardous gases. Consult local regulations and your company's safety procedures.

### Instructions

#### Installation for quick connect fittings

1. Check the package contents for damage. Contact your supplier if any items are damaged or missing.
2. The SPure filter is delivered in the active state, filled with helium. It is ready for use, but it is recommended that the trap be purged with the gas to be purified.
3. Install the quick connect tube fitting ends to the gas lines. For the best connections, start with freshly-cut tubing in the gas line. 1/8" nuts should be turned 3/4 turn past finger tight. 1/4" nuts should be turned 1-1/4 turns past finger tight.
4. Attach the outlet end of the filter by screwing the quick connect fitting nut onto the adapter. This will open both check valves on the outlet side.
5. Attach the inlet end of the filter by screwing the quick connect fitting nut onto the adapter. This will open both valves on the inlet side.
6. Purge the entire system until a total of 1 to 2 liters of gas has been flushed through.
7. Leak check all connections.

## Specifications

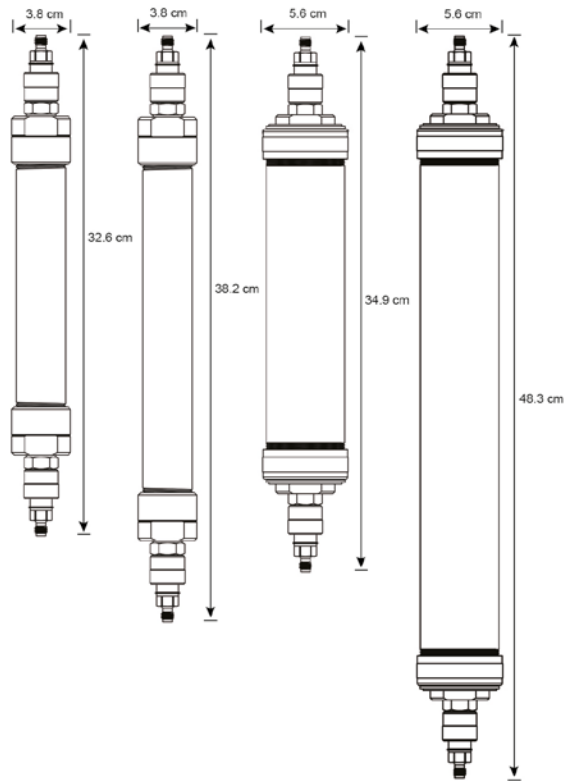


Figure 1. From left to right: 88 cc, 116 cc, 240 cc, 400 cc

## Product Numbers

88 cc	
1/8" Brass	1/4" Brass
202324-B-QC	202325-B-QC
1/8" Stainless steel	1/4" Stainless steel
202324-SS-QC	202325-SS-QC
116 cc	
1/8" Brass	1/4" Brass
202324L-B-QC	202325L-B-QC
1/8" Stainless steel	1/4" Stainless steel
202324L-SS-QC	202325L-SS-QC
240 cc	
1/8" Brass	1/4" Brass
202324D-B-QC	202325D-B-QC
1/8" Stainless steel	1/4" Stainless steel
202324D-SS-QC	202325D-SS-QC

400 cc	
1/8" Brass	1/4" Brass
202324XL-B-QC	202325XL-B-QC
1/8" Stainless steel	1/4" Stainless steel
202324XL-SS-QC	202325XL-SS-QC
Replacement cartridge	
130 cc	
Brass	Stainless steel
202324RP-B-QC	202324RP-SS-QC
475 cc	
Brass	Stainless steel
202324LRP-B-QC	202324LRP-SS-QC
500 cc	
Brass	Stainless steel
202324DRP-B-QCB	202324DRP-SS-QC
750 cc	
Brass	Stainless steel
202324XLRP-B-QC	202324XLRP-SS-QC

## References



### Recycling

Please contact Trajan for information regarding recycling purifier products.

## Information and support

Visit [www.trajanscimed.com](http://www.trajanscimed.com) or contact [techsupport@trajanscimed.com](mailto:techsupport@trajanscimed.com)

Specifications are subject to change without notice.

US AGENT	<b>Trajan Scientific Americas Inc</b>
	2601 Technology Drive,
	Louisville, KY 40299, USA
	Tel: +1 502 491 6300