

ZPure™ Glass H₂O filters

Quick connect fittings

- Do not use without the protective outer plastic tube in place.
- Special precautions might be required when using hydrogen or hazardous gases. Consult local regulations and your company's safety procedures.
- Color Disclaimer: This product is intended to be yellow. Slight variations in shade may occur due to the manufacturing process and differences in individual color perception.

Introduction

Pure gas is critical requirement in gas chromatography, spectroscopy, optics, lithography, and numerous other applications in manufacturing and analytical laboratories. The ZPure™ H₂O filters remove H₂O to low ppb levels.



Important



Warnings must be read carefully and understood. Improper use of this product can cause harm or death to personnel and damage to property!

- Wear eye protection and use caution when working with pressurized systems.
- Maximum pressure is 11 bar / 160 psi.
- When using a high-pressure gas source over-pressure protection must be provided.
- Do not open filter, even after use.
- The rapid oxidation of the active adsorbent in this filter can cause the filter bed to become extremely hot, resulting in danger of fire or burn injuries. Limited exposure of the filter to air during installation does not result in rapid reaction or dangerous heat generation.
- Do not pass air or oxidizing gases in concentrations above 0.1% through this filter.
- Do not expose the protective outer plastic tube to solvents.

Instructions

Installation for quick connect fittings

1. Check the package contents for damage. Contact your supplier if any items are damaged or missing.
2. The filters are delivered in the active state, filled with helium. It is ready for use, but it is recommended that the trap be purged with the gas to be purified.
3. Install the quick connect tube fittings ends to the gas lines. For the best connections, start with freshly-cut tubing in the gas line. 1/8" nuts should be turned 3/4 turn past finger tight. 1/4" nuts should be turned 1-1/4 turns past finger tight.
4. Attach the outlet end of the filter by screwing the quick connect fitting nut onto the adapter. This will open both check valves on the outlet side.
5. Attach the inlet end of the filter by screwing the quick connect fitting nut onto the adapter. This will open the check valves on the inlet side.
6. While maintaining a low purge rate of 10 to 30 cc/min, purge the entire system until a total of 1 to 2 liters of gas has been flushed through.
7. Leak check all connections.

Specifications

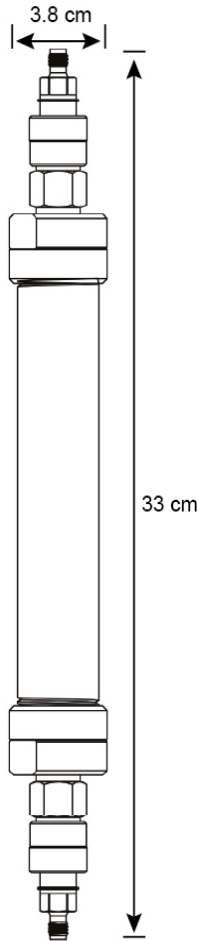


Figure 1. 46 cc

References



Recycling

Please contact Trajan for information regarding recycling purifier products.

Information and support

Visit www.trajanscimed.com or contact techsupport@trajanscimed.com

Specifications are subject to change without notice.

US
AGENT

Trajan Scientific Americas Inc
2601 Technology Drive,
Louisville, KY 40299, USA
Tel: +1 502 491 6300

Product numbers

130 cc	
1/8" Brass	1/4" Brass
202234-B-QC	202235-B-QC
1/8" Stainless steel	1/4" Stainless steel
202234-SS-QC	202235-SS-QC
Replacement cartridge	
130 cc	
Brass	Stainless steel
202234RP-B-QC	202234RP-SS-QC