

Introduction

Pure gas is critical requirement in gas chromatography, spectroscopy, optics, lithography, and numerous other applications in manufacturing and analytical laboratories. The ZPure™ H₂O filters remove H₂O from inert gases, He, Ar, N₂, H₂ methane and clean dry air (CDA) to low ppb levels with very high capacity.



Important



Warnings must be read carefully and understood. Improper use of this product can cause harm or death to personnel and damage to property!

- Wear eye protection and use caution when working with pressurized systems.
- Maximum pressure is 13.8 bar / 200 psi.
- When using a high-pressure gas source over-pressure protection must be provided.
- Do not open filter, even after use.
- Do not use with supply gas H₂O concentration greater than 0.1%. Contains molecular sieve which may become very hot in contact with liquid water. Do not use with liquids.
- Special precautions might be required when using hydrogen or hazardous gases. Consult local regulations and your company's safety procedures.

Instructions

Installation for quick connect fittings

1. Check the package contents for damage. Contact your supplier if any items are damaged or missing.
2. The filters are delivered in the active state, filled with helium. It is ready for use, but it is recommended that the trap be purged with the gas to be purified.
3. Install the quick connect tube fittings ends to the gas lines. For the best connections, start with freshly-cut tubing in the gas line. 1/8" nuts should be turned 3/4 turn past finger tight. 1/4" nuts should be turned 1-1/4 turns past finger tight.
4. Attach the outlet end of the filter by screwing the quick connect fitting nut onto the adapter. This will open both check valves on the outlet side.
5. Attach the inlet end of the filter by screwing the quick connect fitting nut onto the adapter. This will open the check valves in both the quick connect tube fitting and the quick connect adapter, allowing gas to flow through.
6. Purge the entire system until a total of 1 to 2 liters of gas has been flushed through.
7. Leak check all connections.

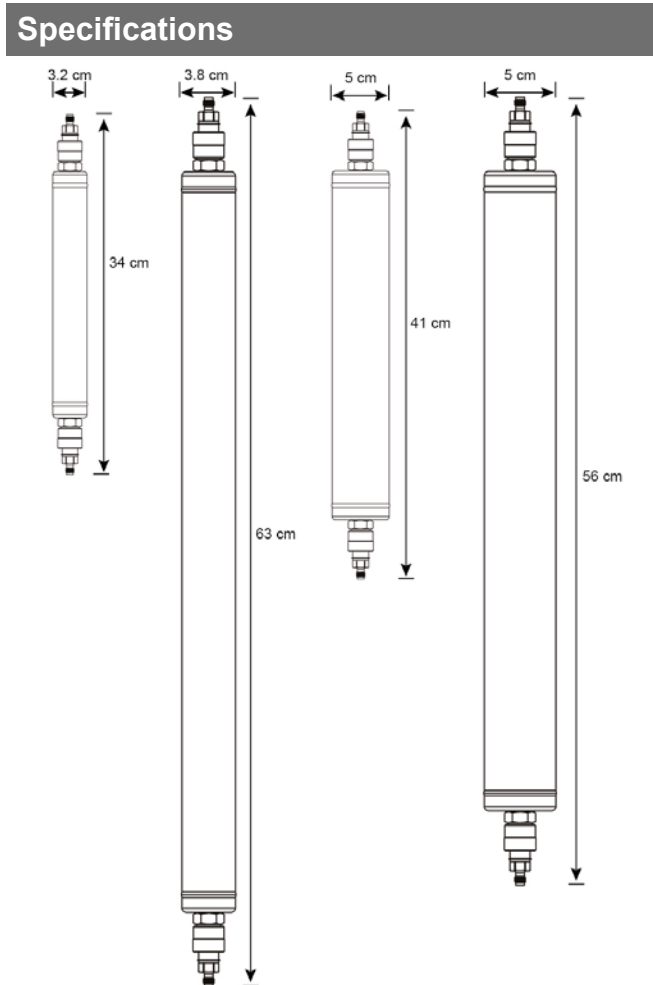


Figure 1. From left to right: 130 cc, 475 cc, 500 cc, 750 cc

Product numbers

130 cc	
1/8" Brass	1/4" Brass
202230-B-QC	202231-B-QC
1/8" Stainless steel	1/4" Stainless steel
202230-SS-QC	202230-SS-QC
475 cc	
1/8" Brass	1/4" Brass
202230L-B-QC	202231L-B-QC
1/8" Stainless steel	1/4" Stainless steel
202230L-SS-QC	202231L-SS-QC
500 cc	
1/8" Brass	1/4" Brass
202230D-B-QC	202231D-B-QC
1/8" Stainless steel	1/4" Stainless steel
202230D-SS-QC	202231D-SS-QC

750 cc	
1/8" Brass	1/4" Brass
202230XL-B-QC	202231XL-B-QC
1/8" Stainless steel	1/4" Stainless steel
202230XL-SS-QC	202231XL-SS-QC
Replacement cartridge	
130 cc	
Brass	Stainless steel
202RH2-B-QC	202RH2-SS-QC
475 cc	
Brass	Stainless steel
202RH2L-B-QC	202RH2L-SS-QC
500 cc	
Brass	Stainless steel
202RH2D-B-QCB	202RH2D-SS-QC
750 cc	
Brass	Stainless steel
202RH2XL-B-QC	202RH2XL-SS-QC

References



Recycling

Please contact Trajan for information regarding recycling purifier products.

Information and support

Visit www.trajanscimed.com or contact techsupport@trajanscimed.com

Specifications are subject to change without notice.

US AGENT	Trajan Scientific Americas Inc
	2601 Technology Drive,
	Louisville, KY 40299, USA
	Tel: +1 502 491 6300